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Guangzhou Huayuan Pharmaceutical Equipment Co., Ltd

地址：广东广州市番禺区大石街道石南1路潮联创业中心225
No.225 ChaoLian Venture Center, Shinan 1 road, Dashi street,
Panyu District, Guangzhou City, Guangdong Province

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“tandardized management, safe Huayuan”

safety production construction policy;

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Huayuan people's goal is to

"survive with quality and drive by science and technology";

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The service tenet of

"receiving feedback in good faith and serving warmly and actively in time";

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Understand the new value of enterprises,

has been for the development of regional economy

and people's health and strive!

About Us

Guangzhou Huayuan Pharmaceutical Equipment Co., Ltd. is an innovative technology enterprise specializing in pharmaceutical equipment, biological fermentation equipment, plant extraction equipment, vacuum low-temperature drying equipment.

The company's existing mature products: multi-functional extraction tank, extraction and concentration equipment, evaporation and concentration equipment, alcohol recovery equipment, filtering equipment, crystallization equipment, container equipment, pump valve fluid pipe fittings and other products.

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GuangDong Guangzhou Huayuan Pharmaceutical Co. LTD. a special manufacturer integrating design, development, manufacture, engineering construction and after-sale service, devotes itself to manufacture and sale the complete equipment for pharmacy

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The company owns a team of excellent specialists in design and manufacture. It mainly produces health-grade equipment for extraction, concentration and reclamation and stainless steel tank and pipe accessories and valve, series full automatic CIP cleaning machine and PLC system. The product is designed and made strictly in accordance with the ISO90012000, meeting with the GMP standard.

The union, enterprise and high talent is the base of our success. Our advanced craft in extraction and concentration of China traditional drug and reclamation of alcohol and high efficient and energy-saving equipment and the high practice and attractive

appearance can improve the production efficiency greatly; The complete equipment for large transfusion and water injection features pleasant appearance and special structure;The injection and material allocation system and extraction & concentration system for China traditional drug and the CIP system can realize the automatic management for production and control.

The HuaYuan company holds the marketing purpose of WIN-WIN and keeps long-term cooperation with the famous manufacturer home and abroad. The product is exported to Europ,Australia and Southeast Asia and enjoys high appreciation from the client.

因为 华远 始终坚持“真诚接受反馈意见、主动及时热情服务”的服务宗旨 实现“华远-客户”双赢。
Because Huayuan always adhere to the "sincere acceptance of feedback, proactive and timely warm service" service purpose, to achieve the "Huayuan - customer" win-win

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Accord with FDA standard, meeting with the GMP requirement, ISO9001 certification, first-class and second-class pressure vessel manufacture.

Perfect quality originates from advanced manufacturing methods

High-technology and advanced skills for guarantee the excellent quality and performance.

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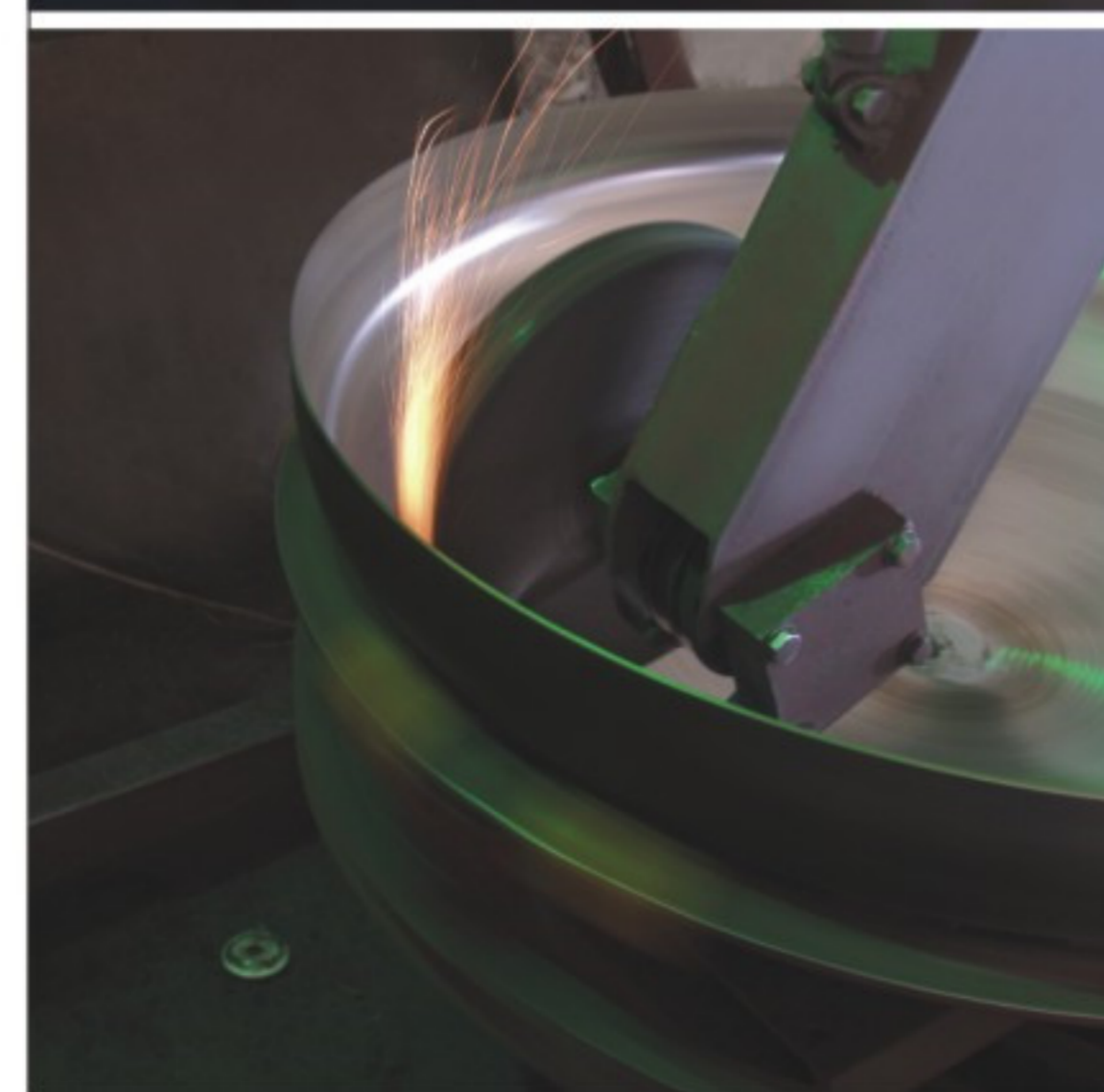
Our factory has equipped advanced TIG welding machine, various manual and automatic welding machine, large-scale closure head polishing set, spinning machine and bending machine, CNC etc. we also offers non-standard equipment in special design that can match your highly demand.

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The success of HuaYuan is not only the perfect of the surface, but also the fineness of details. We strictly examine every fitting, every craft, and every machine, and turn them into the representative of the finest once it is taken out of our factory.

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Advanced automatic production equipments, steam line production, matched first-class quality. Chasing the perfection is our spirit.

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"Unstoppable pursuit of high quality for the customers" that are HuaYuan's main line, as a well-known machinery manufacturers, we believe that only chasing perfect quality is the constant development of the enterprise. The company has modern production and testing equipments and strictly following the ISO9001 Quality Management System Requirement. "Excellent performance and perfect quality is the operating guidelines of HuaYuan. We will produce first-class products, make perfect service for customers!"

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Project Examples

Pharmaceutical Equipment

We have developed pharmaceutical machinery and equipment that can not only meet your requirement of safe, efficient, and sanitary, But also ensure the detailed processing of the product.

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Herb extraction concentration complete sets system; infusion, water injection system of complete sets equipment; biological fermentation complete sets equipment, suger processing system, raw materials, equipment Synthetic Medicine; Pet medicine products equipment; emulsifying system equipment;

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TQ series multi-functional Extractor

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- ◆Application: This device is applied in industries such as herb, food and chemistry etc. Such as normal pressure, low pressure, water decocting, immersion, thermal circumfluence, compelled recycle and lixiviation, extraction of aroma oil etc.
- ◆Characteristics: High efficiency and convenient operation
- ◆Configure: CIP cleaning system
- ◆Standard: Matched with GMP

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
- ◆Security design: The latest residue outlet with high security, it open with pneumatic cylinder and lock with dual-gas cylinder accompanied with a secure lock cylinder. The complete device is easy to open and operate, and it has a reliable self-lock performance.
- ◆This control eliminates the leakage or detachment caused from the unstable gas pressure different with the others

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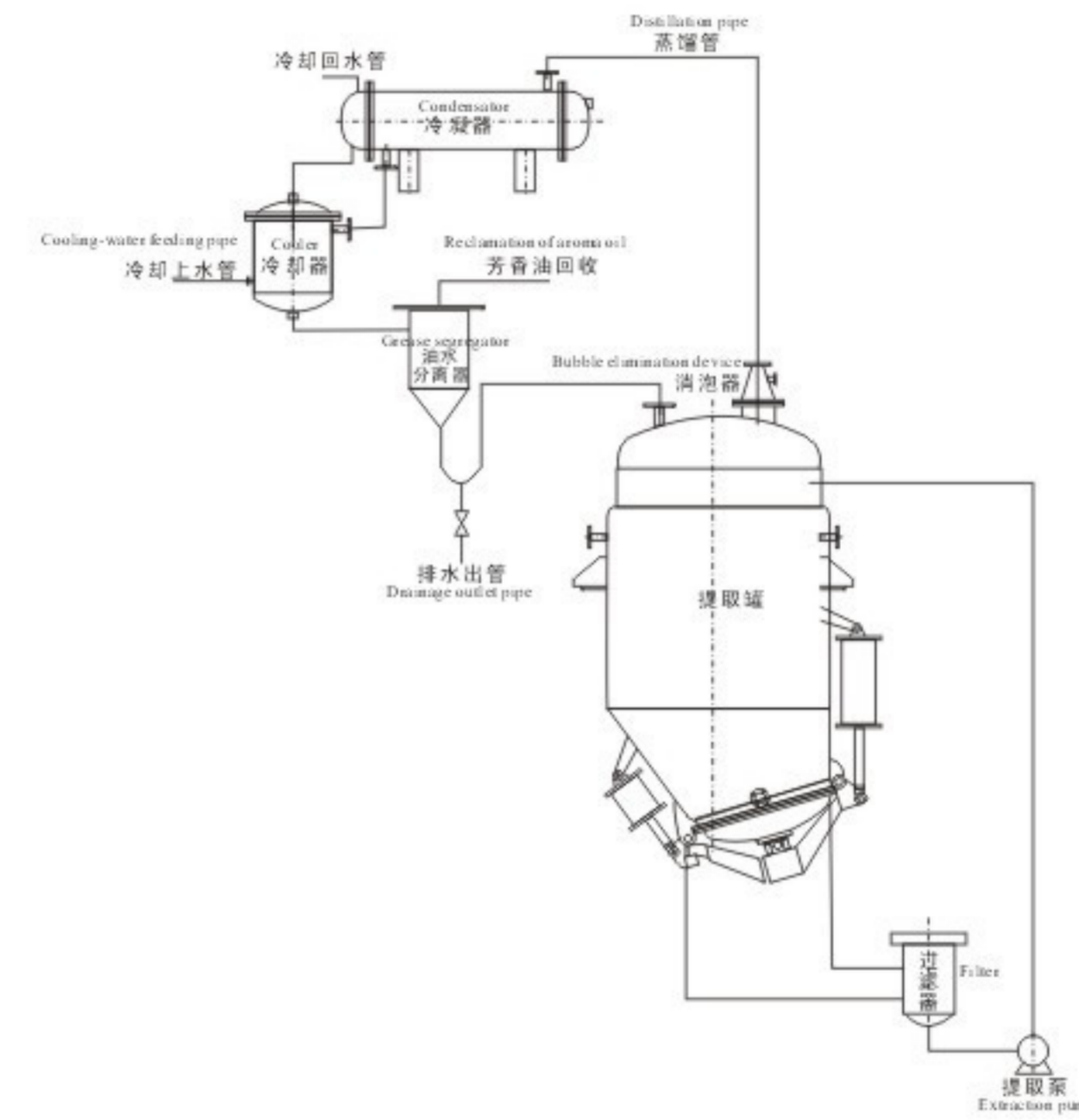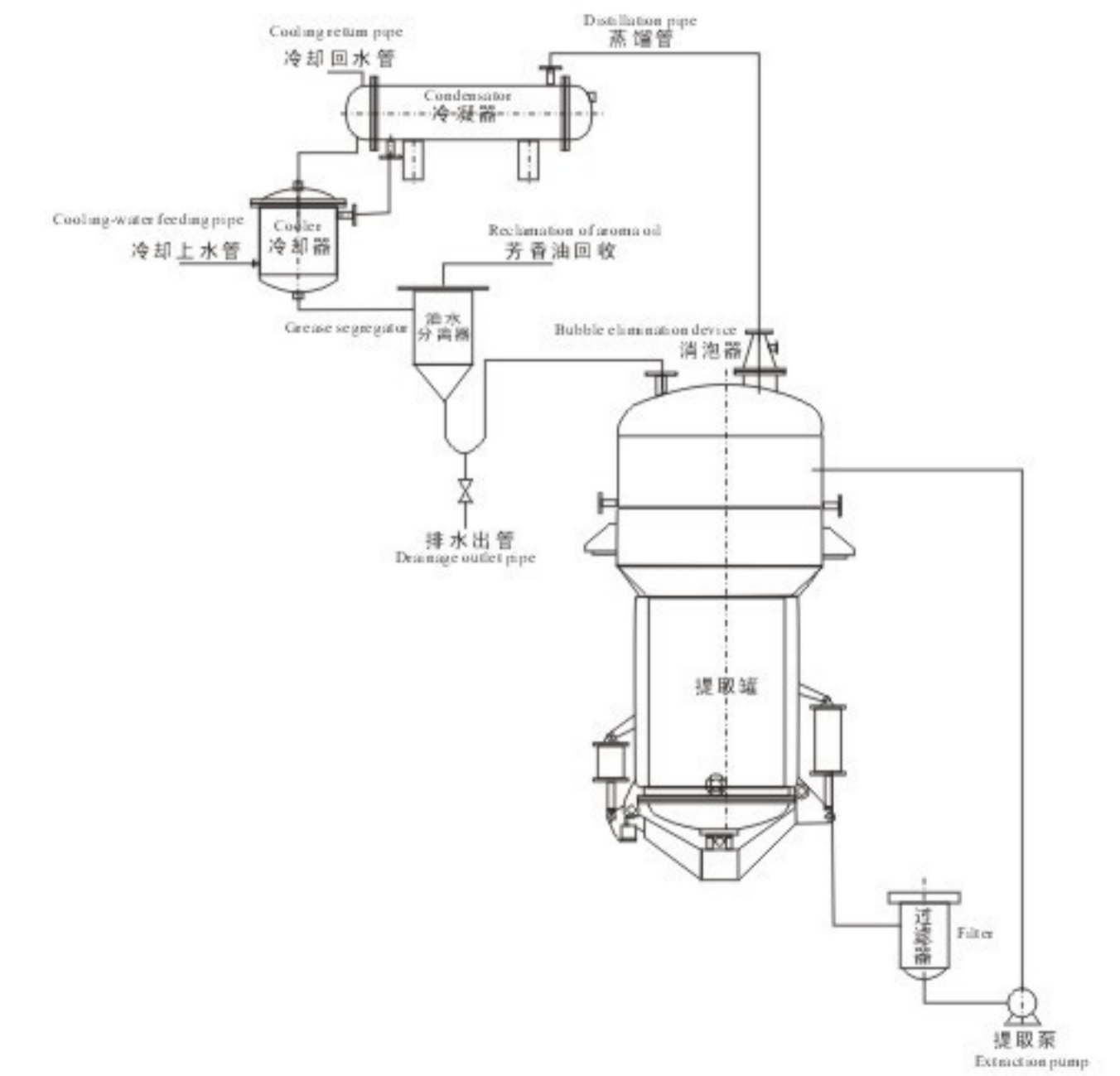
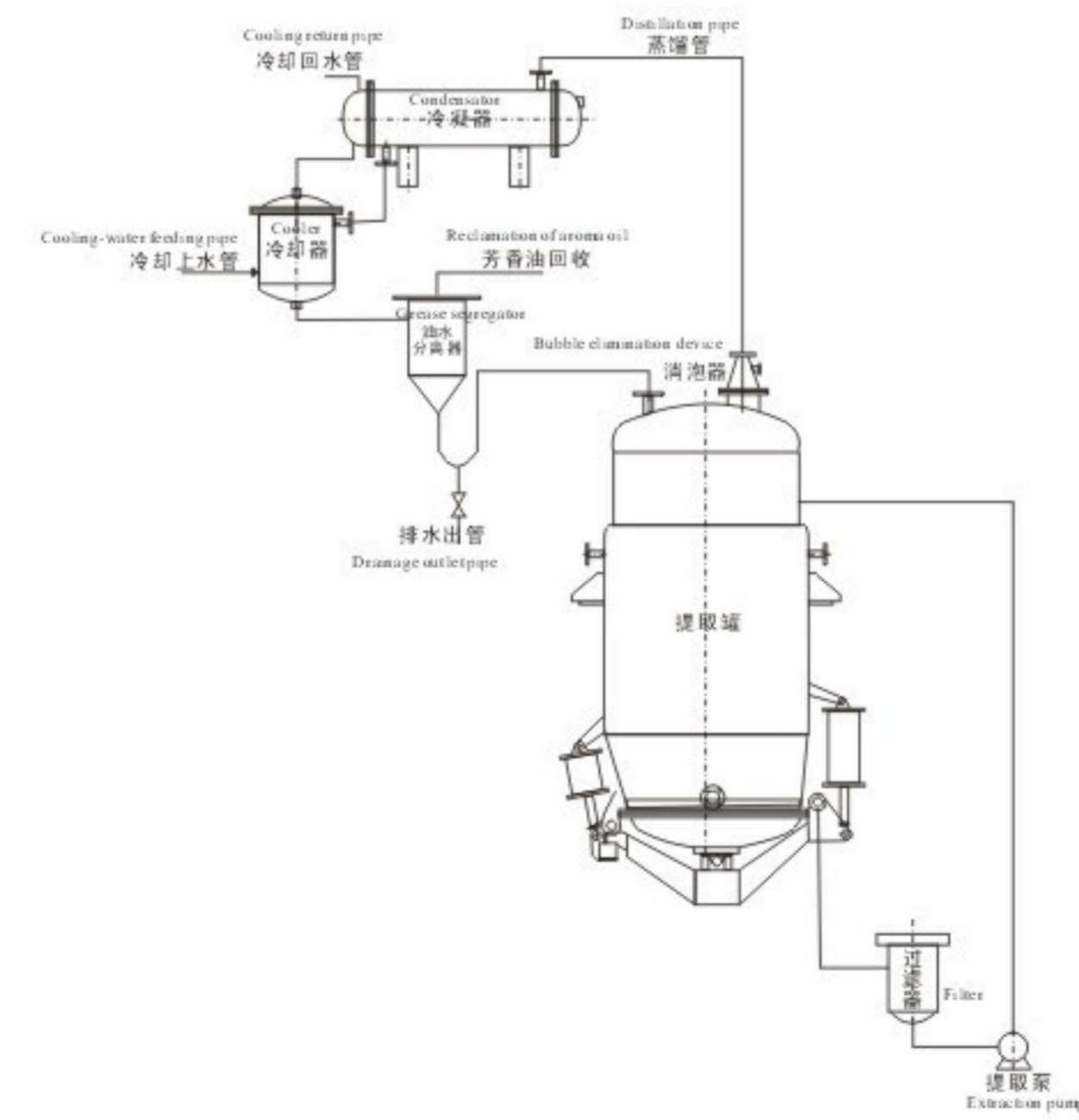
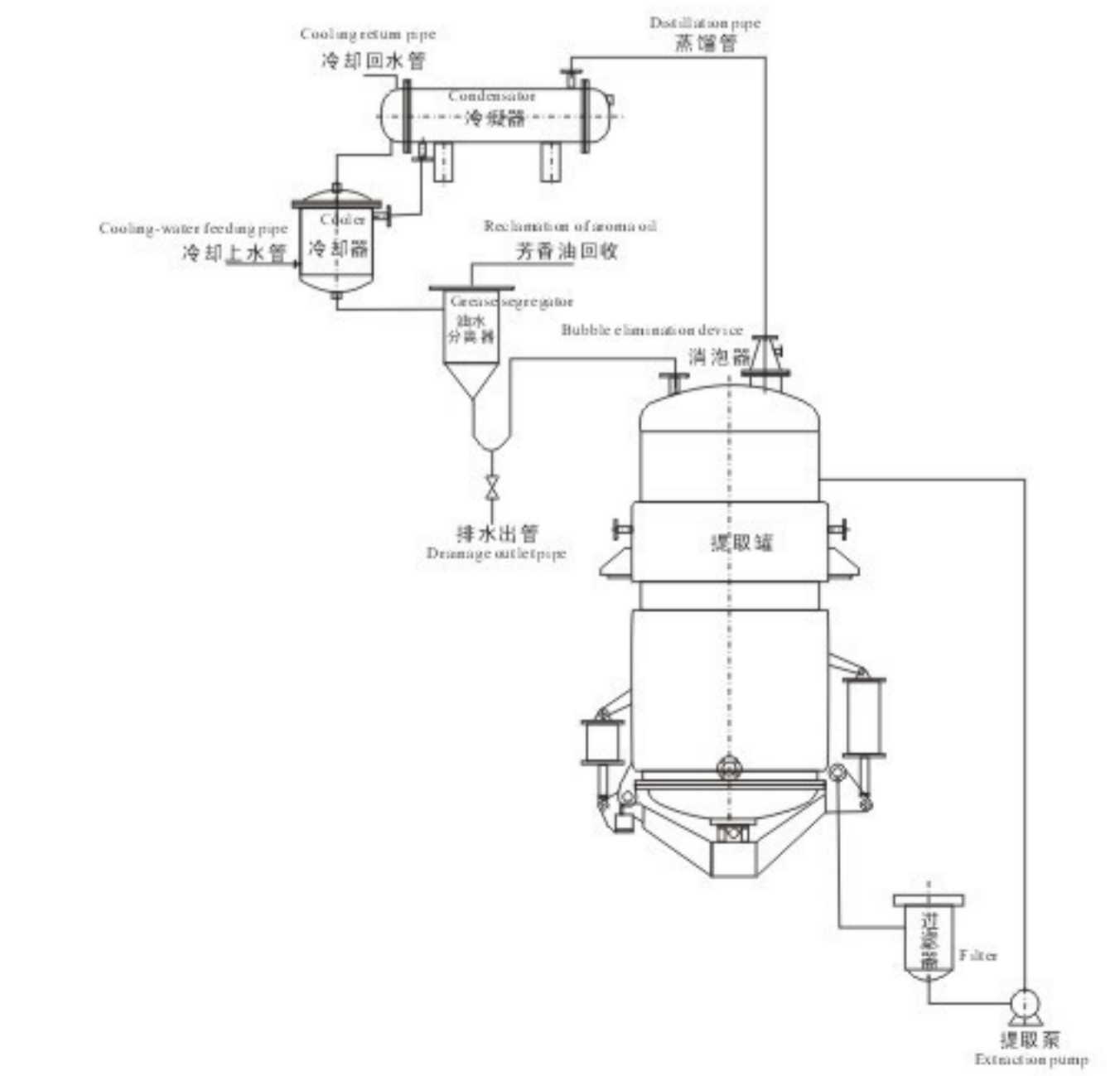
DTQ series multifunctional extractor (dynamic agitator)

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TQ-D 直锥型(带搅拌)
TQ-D Vertical Prick Type(with milling)

- ◆Application: This device is applied in industries such as herb, food and chemistry etc Such as normal pressure, low pressure, water decocting, immersion, thermal circumfluence, compelled recycle and lixiviation, extraction of aroma oil etc
- ◆Spiral agitator and propel agitator
- ◆Characteristics: High efficiency and convenient operation
- ◆Configure: CIP cleaning system
- ◆Standard: Matched with GMP

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- ◆Energy-saving set
- ◆Integrated with 3 functions that simplified the operation procedure.
- ◆under the close and seal condition, no leakage
- ◆SUS316L(or SUS304), meeting with the GMP standard.
- ◆best applied in hospital, university laboratory, pharmaceutical factory and R&D institute and food manufacture to extract and concentrate the herb and juice etc

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The device applied in herbal medicine, China traditional drug, pet drug, food additive etc, to extract with water, ethanol, carbinol and acet.

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Equipment Characteristics

◆ The rate for cream collection is 10~15% higher than that of the multifunctional tank, and the effective content inside the cream is one time higher than the normal one. Because the heat solvent is added into the drug surface continuously and then they pass the drug layer from up to down with high speed to solve the solute in the drug material during the extraction, so the solute in the drug material and the solute in the solvent keep high grads. If the solute in the drug material can separate out with high speed until it separates out, then the extraction extent of the effective content is high. Because the concentration is finished inside the seal device with little loss and high transformation rate, so the rate for cream collection is high and the effective content inside the drug cream is high.

◆ For the factors of high speed of separation out of solute, short time of extraction synchronous operation of concentration and extraction, the time from changing the drug material into the concentrated drug cream is only seven to eight hours, which highlights high equipment utility.

◆ The equipment is simple and has little use area and features high equipment utility, which highlights economic investment.

◆ The extraction just needs one time of use of solvent, which is used cyclically in a set of seal equipment. The solvent in the drug residue can be reclaimed basically, which makes the solvent quantity is 30% less than that of the multiple function extraction tank and the solvent consumption is 50~70% less than that of the multiple function extraction tank. The above advantages make it available the extraction of organic solvent and purification of effective content in the China traditional drug.

◆ For the concentrated secondary vapor serving as heat source, the extraction liquid into the concentrator is at the same temperature level with that of the concentrator, which can save the vapor for more than 50% and feature easy operation and reduces the operator.

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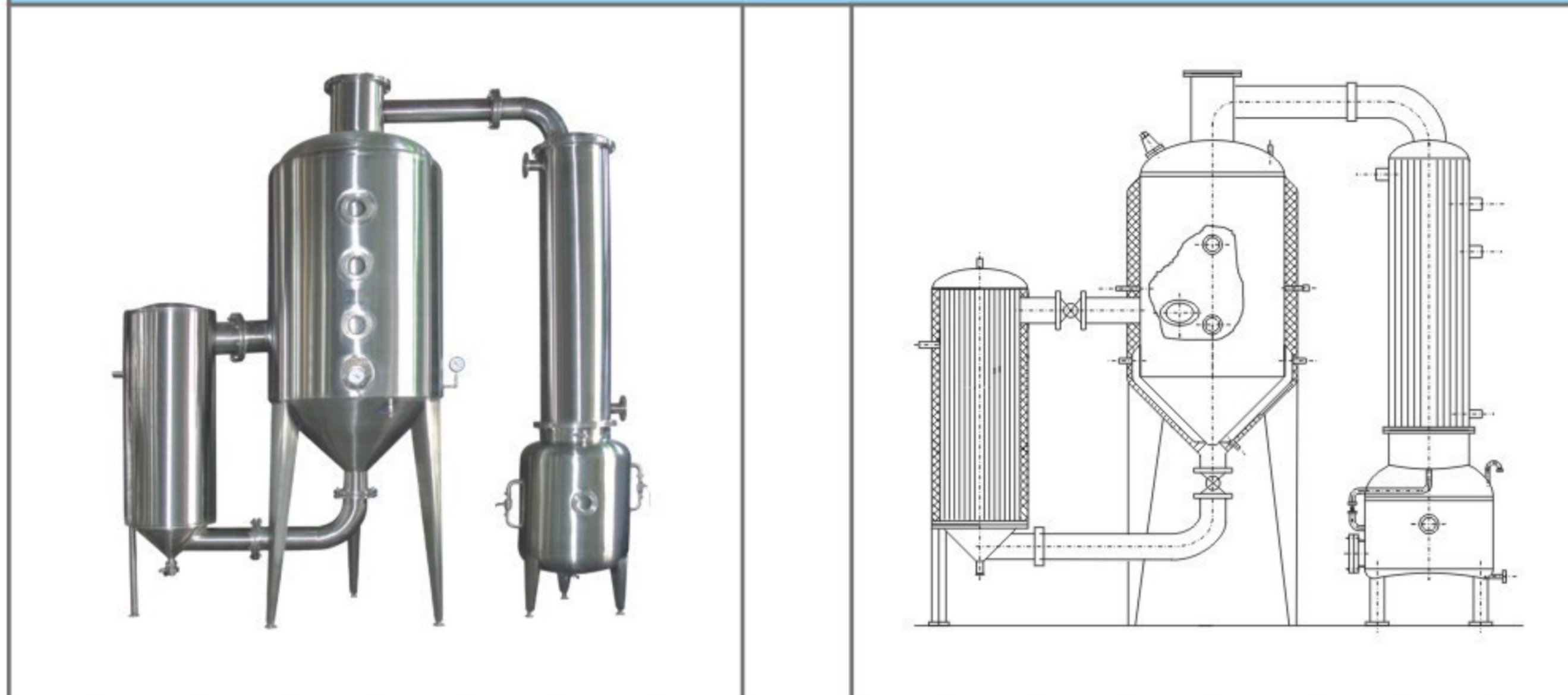
Principle

Put the drug material into the extraction tank and insert the solvent such as water, ethanol, carbinol and acet, etc (in accordance with the technology requirement) whose capacity is five to ten times the drug material. The tee in the extraction tank and the vapor valve of clip sleeve open to heat the extraction liquid until boil that will last for about twenty to thirty minutes. And then, the extraction & filter pipe sorbs one third extraction liquid into the concentrator. And the tee in the extraction tank and the vapour valve of clip sleeve are closed and the heater valve is opened to concentrate the material liquid. The second vapor produced during concentration can enter the extraction tank through the upward pipe at the evaporator to serve as the heat source for extraction to maintain the boil inside the tank.

The second vapor continues rising to change into the condensation liquid by the condenser and falls back inside the extraction tank to serve as new solvent to be added at the drug surface. The new solvent arrives down at the bottom of the extraction tank throughout the drug material layer. The dissolvable effective content at the drug material is resolved into the solvent inside the extraction tank. The extraction liquid enters the concentrator by the extraction & filter pipe and the second vapour produced for concentration arrives the extraction tank to serve as heat source and new solvent. The extraction of new solvent with large circumfluence can make the density of the drug material and solvent at high grads. The solvent in the drug material separates out at high speed until a complete separation (i.e. the extraction liquid is colorless.). Then, the extraction liquid stops entrance of the concentrator. The concentrated second vapor enters the cooler to continue concentration until the drug cream with required proportion (1.2~1.34) is available. The colorless liquid inside the extraction tank can be stored at the storage tank for reuse. The drug residue is drained out of the residue outlet. In the case of extraction with organic solvent, it is necessary to add proper water and open the tee and the clip sleeve vapor valve. After reclamation of solvent, the residue is drained.

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The concentrator is suitable to concentrate the China traditional drug, west drug, dextrose, amylum, monosodium glutamate, food and milk product and chemical material, especially suitable to concentrate the thermal sensitive material under low temperature and vacuum. After our improvement, the feature and profit is much more obvious. The client can choose the various concentrator in accordance with the concentration content.

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Feature & profit

- ◆ Energy-saving profit: In case of SJN-1000, the saves respectively in the terms of vapor, water and electric per year are about 3500 tons, 90000 tons and 80000 watt, equal to RMB 100,000-150,000 Yuan.
- ◆ The device uses the modes of outer-heated natural cycle and vacuum negative pressure evaporation featuring swift evaporation with concentration proportion of 1.2-1.35 (in case of normal China traditional immerse cream) without need of open-mouth pan for collection of cream.
- ◆ The concentrator uses dual-effect to synchronously vapor, which to utmost extent utilize the second vapor. The client can lower investment in boiler and the energy consumption can be 50% less than that of the single-effect concentrator. It is about less than one year that the client can withdraw the investment in the concentrator.
- ◆ Feature of multiple function handle: (1)The reclaimable alcohol concentration is about 80%. (2)The single-effect and dual-effect can time after time collects cream together. (3)The intermittent and continuous material feeding is available.
- ◆ The special structure of concentrator can bring the concentrated material thick taste and even cream quality. The concentration under the seal and non-bubble condition can prevent material loss and pollution. The present material can be concentrated presently, which features immune to block and convenient cleaning which can be make just by opening the up and down covers and can be make by one person.
- ◆ The part inside the concentrator in contact with the material is made with stainless steel SUS316L (or SUS304). The thermal layer made with polyurethane is available at the heater and evaporator. The surface is treated with sand grinding and surface polishing, meeting with the GMP standard.

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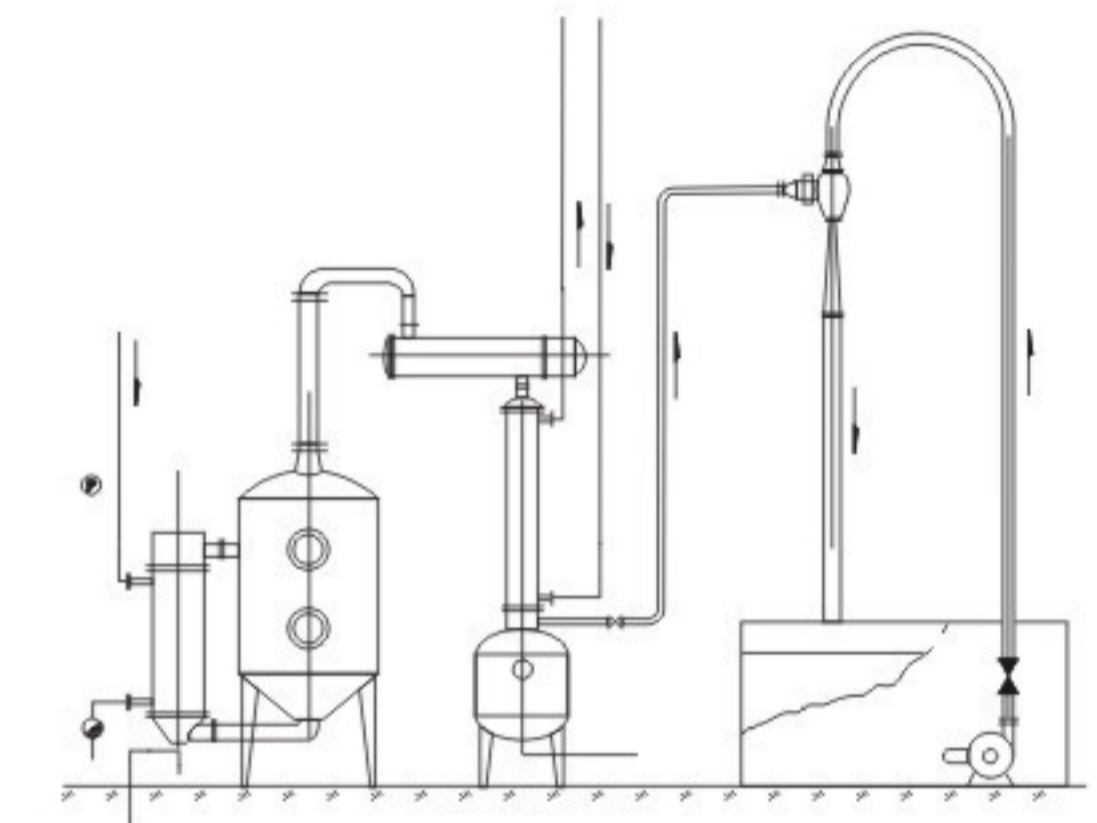
This concentrator is suitable for concentrating the fluid materials of traditional medicine, western medicine, grape sugar, starch, gourmet powder, food and milk products, chemical industry, especially for thermal sensitive matter low temperature vacuum concentration. Our factory has carried out the careful accumulation and meticulous design and reform and got more profits than the original characteristic. You can select the series concentrator in technical parameters according to your concentration degree.

Feature & profit

- ◆Energy-saving profit: Take SJN-1000 as example, the saves respectively in the terms of vapor, water and electric per year are about 5400 tons, 140000 tons and 130000 watt, equal to RMB 200000 Yuan around.
- ◆The device uses the modes of outer-heated natural cycle and vacuum negative pressure evaporation featuring swift evaporation with concentration proportion of 1.4 (in case of normal China traditional immerse cream).
- ◆The concentrator uses triple-effect to synchronously vapor, which to utmost extent utilize the second vapor. The client can lower investment in boiler and the energy consumption can be 70% less than that of the single-effect concentrator. It is about less than one year that the client can withdraw the investment in the concentrator.
- ◆Feature of multiple function handle: 1 The reclaimable alcohol concentration is about 80%. 2 The triple-effect can concentrate three different materials. 3 The single-effect, dual-effect and triple-effect can collect cream together time after time. 4 The intermittent and continuous material feeding is available. 5 The one-off cream formation is available. The material input is finished by the first effect and after the second and third effect, the finished cream is available one-off.
- ◆The special structure of concentrator can bring the concentrated material thick taste and even cream quality. The concentration under the seal and non-bubble condition can prevent material loss and pollution. The present material can be concentrated presently, which features immune to block and convenient cleaning which can be make just by opening the up and down covers and can be make by one person.
- ◆The part inside the concentrator in contact with the material is made with stainless steel SUS316L (or SUS304). The thermal layer made with polyurethane is available at the heater and evaporator. The surface is treated with sand grinding and surface polishing, meeting with the GMP standard.

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JN multi-functional alcohol reclamation concentrator

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◆Alcohol reclamation: The large reclamation capacity and the vacuum operation mode improve the craft process with a reclamation rate of alcohol of 95% and a loss rate of alcohol below 5%. The power consumption lowers for 30%, which improves that the device is a energy-saving equipment with economical investment and high reclamation economy profit.

◆Cocentrated liquid material: The device uses the modes of outer-heated natural cycle and vacuum negative pressure evaporation featuring swift evaporation with concentration proportion of (1.4); The liquid material concentrates without bubble under full seal state. The drug liquid after concentration by the device features non-contamination and thick drug taste and easy cleaning(it is just to open the up and down cover for cleaning.). The device is easy to handle and it needs a little foundation area for installation. The heater and Evaporator use stainless steel SUS316L (or SUS304) for thermal purpose. The polyurethane or pearl cotton is used as the thermal material. The surface is treated with sand grinding and surface polishing, meeting with the GMP standard.

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
The device is to concentrate the material in the occasions such as China traditional drug, west drug, dextrose, amyllum sugar, oral liquid, chemistry, foodstuff, monosodium glutamate and vintage and to reclaim the industrial organic solvent such as alcohol. It is suitable to concentrate the thermal sensitive material under lower temperature and vacuum with little batch number and multiple varieties.

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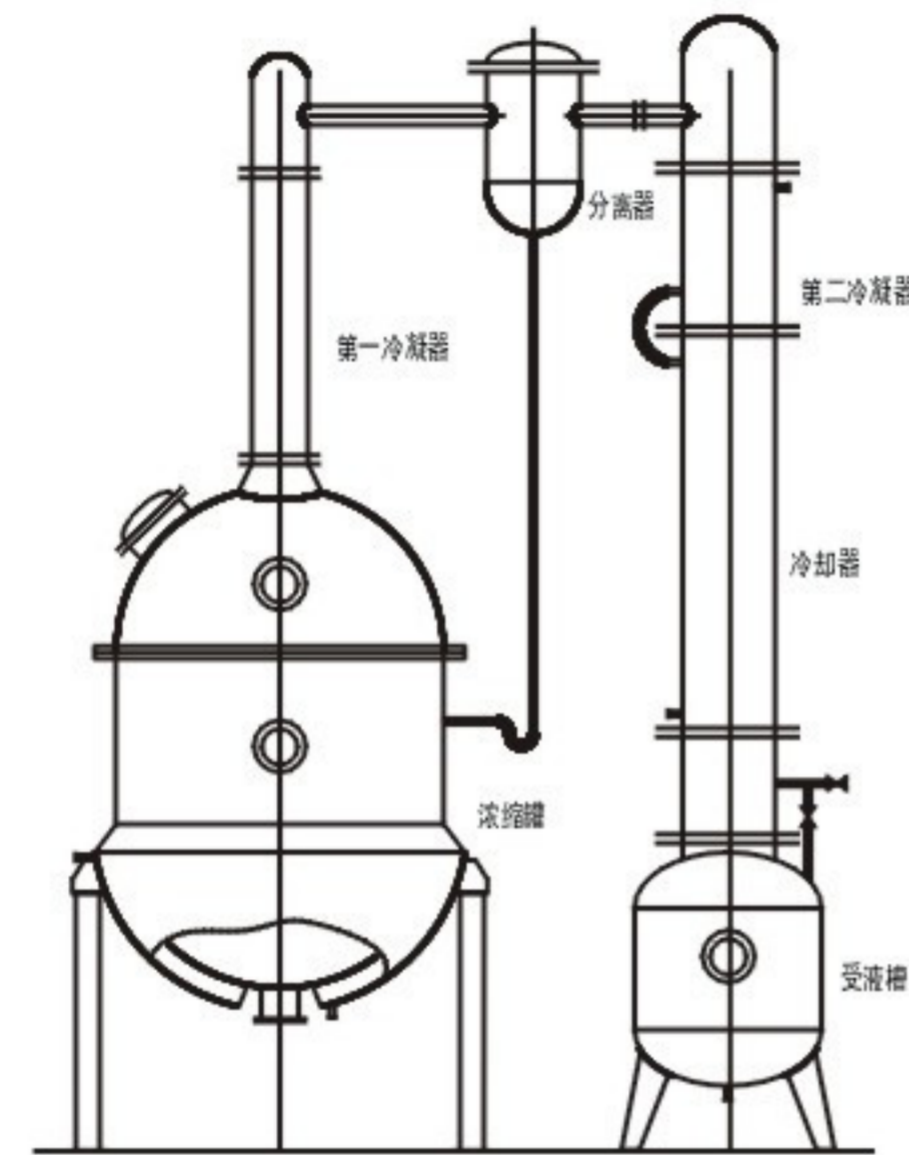ZN vacuum pressure reduction concentrator

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
The device includes concentration tank, condenser, gas-liquid separator, cooler and liquid receiving barrel. The concentrator tank is of clip sleeve structure. The condenser is of line pipe type and the cooler is of tray pipe type. The device is suitable to concentrate in the pharmacy, foodstuff and chemistry and also can serve as reclamation of alcohol and simple reflux extraction. The fact that the part in contact with the material is made with stainless steel SUS316L (or SUS304) brings qualified anti-corrosion performance, meeting with the GMP standard.

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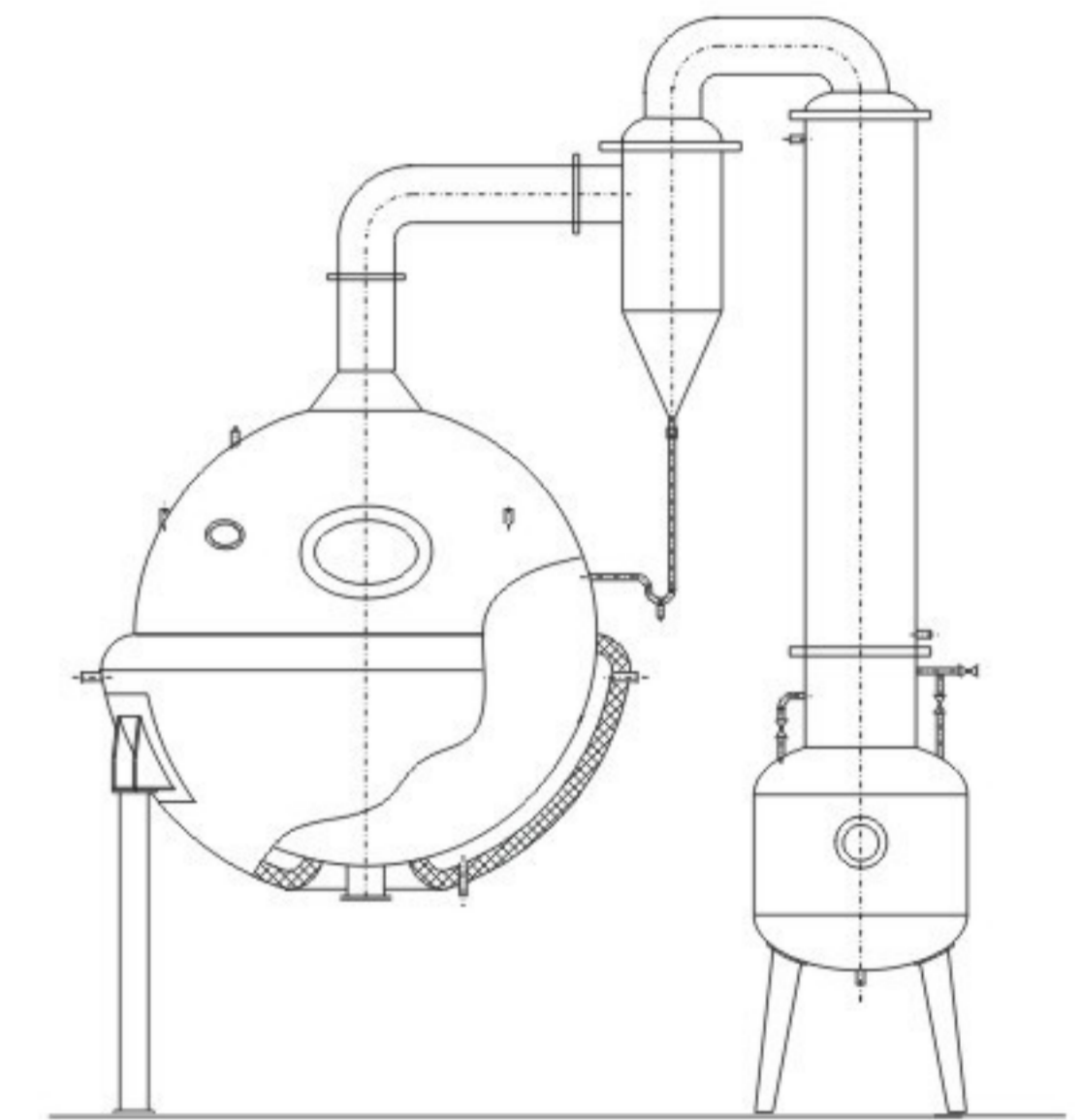The roundness concentration tank, consisting of main body, condenser, gas-liquid separator and liquid receiving barrel, serves to concentrate, distil the material liquid and reclaim organic menstruum in the pharmacy, foodstuff and chemistry. The concentration under reduced pressure brings short time of concentration and prevent the destruction of the effective content in the thermal sensitive material. The fact that the part in contact with the material is made with stainless steel SUS316L (or SUS304) brings qualified anti-corrosion performance, meeting with the GMP standard.

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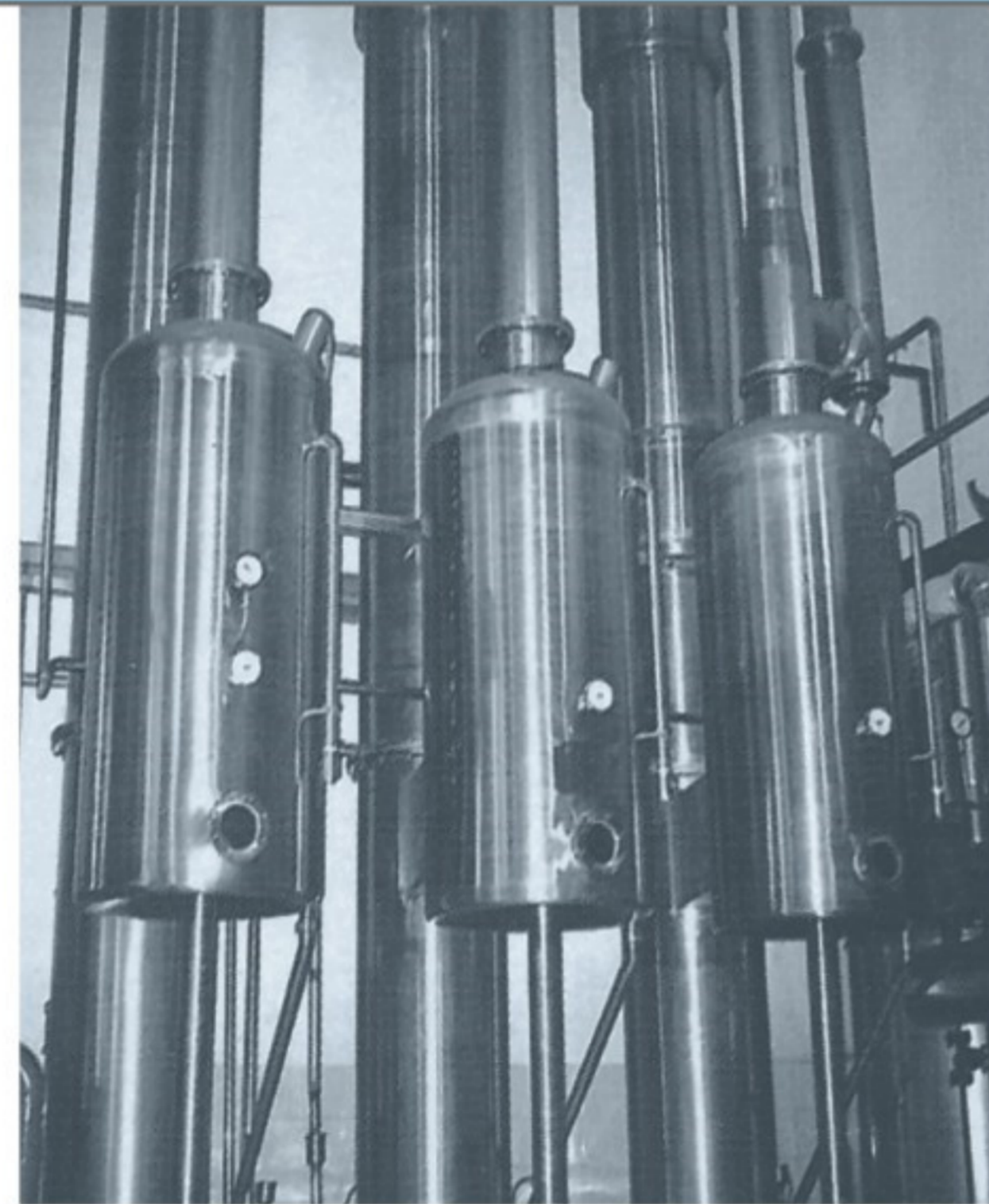
LJN Three effect falling-film continuous feeding concentrator

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
Accord with FDA standard,meeting with the GMP requirement,ISO9001 certification, first-class and second-class pressure vessel manufacture.

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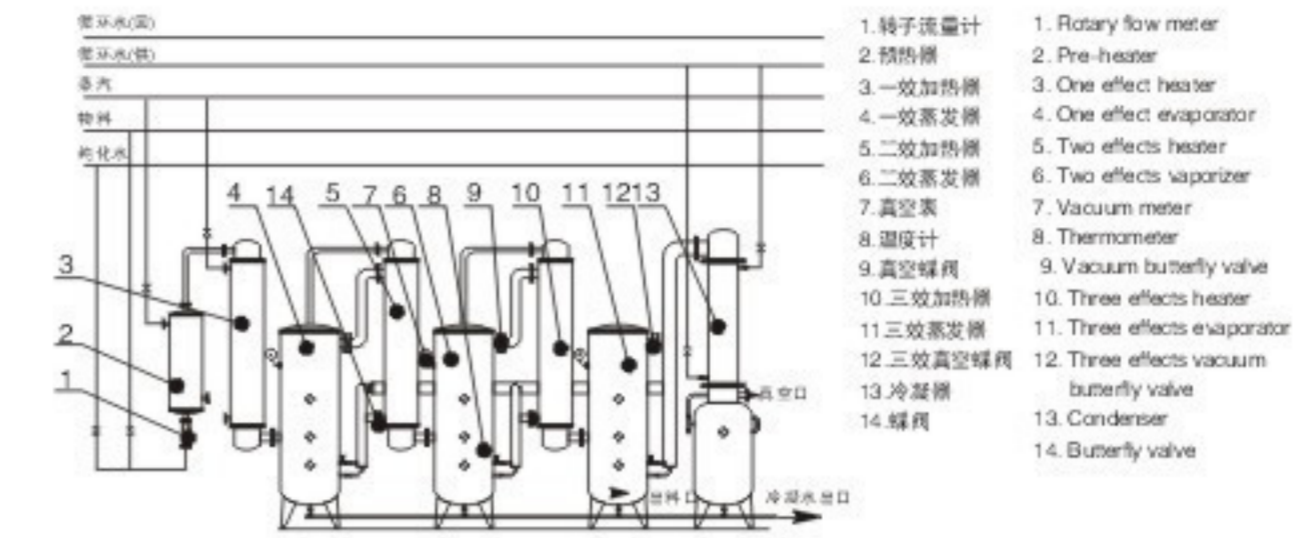The presently used concentrator began serving at the end of 1980's and the Beginning of 1990's, with obvious shortcoming at application:

- ◆ One concentration cycle is of two times of feedings, with different discharge concentration (density) of three effects. The material concentration of three effects is usually lower. So it is a must to mix first effect concentration or first effect concentrated material to meet technology requirement, but which is difficult to be under control.
- ◆ Complicated operation. The feeding is of intermittent type. It is necessary to send the three effect material to the first effect and second effect, which leads to the density of concentration material difficult to control and is liable to produce wrong operation to cause loss to production.
- ◆ Narrow adaptation to material. The structure mode is outer recirculation flat flow feeding. The material leaves the difference of density as push force for recirculation and obtains density difference recirculation by repeated heating. The coefficient of heat transmission is slightly low and the temperature difference for heating is somewhat large(10-15℃), meanwhile, the small push force to material flow results in slow flow speed. The adaptation scope to material viscosity is narrow (0-100cp), destroy to effective element of heat-sensitive material is rather serious and the material is bound to be crystallized, then to deposit and even coke.
- ◆ Low energy-saving effect. The low heat transmission coefficient leads to low heat efficiency that can't save energy very much. We integrated the highlight of imported concentration equipment with that of the counterparts used at other industry and the characteristic of extraction liquid of Chinese drug to introduce the client series LJN type three effect (double effect, single effect) falling-film continuous feeding concentrator. The equipment feature is as follows:
- ◆ Continuous feeding and discharge, stable and convenient operation. The only operation package of setting of different parameters at the beginning of feeding, assurance of reliable and stable feeding supply and easy regulation can certainly produce continuous, stable and efficient concentration, get qualified density and simultaneously recover solvent.

◆ Large concentration density, (1.25-1.28), high heat efficiency and effective energy-saving fruit. To this system, we used optimum to leave a vaporization heat transmission coefficient of 3400-3800 w(m²·K), which is very benefit to concentration of Chinese extraction liquid. The cause for it is that the material flows quickly with the form of film inside the acceleration pipe, and simultaneously, the temperature difference for heat is somewhat small, (just 8-10℃). The material can form quick flashing, the medicine liquid stays and be heated for short time; the adaptation scope of material viscosity is wide, (50-500cp); the even heating excludes crystallization, and then to deposition and even coke. So it features little destroy to effective element, full utilization of heat energy and fruitful energy-saving.

◆ Convenient to employment of advanced control mode. The temperature index for series LJN concentration flow is mainly subject to the temperature of each of effect. Each of parameter keeps stable during system performance, which benefits the utilization of advanced control unit featuring (PLC and man-machine interface control) and leaves the lower cost of control equipment, makes the production of Chinese drug by automatic control great convenience.


◆ Advanced and reasonable structure. We used optimum design scheme. Advanced distribution device is responsible for liquid is attribution during feeding and flow inside the pipe to make the material able to enter heating tube evenly. Simultaneously, the heating tube, made by qualified and imported inner-wall mirror tube, can reduce the flow resistance to the material inside the tube, increase the flow speed, and prevent crystallization and deposition. To lift concentration efficiency, the advanced mist collector can eliminate mist. The separation efficiency can be of 96-99%, with low resistance.

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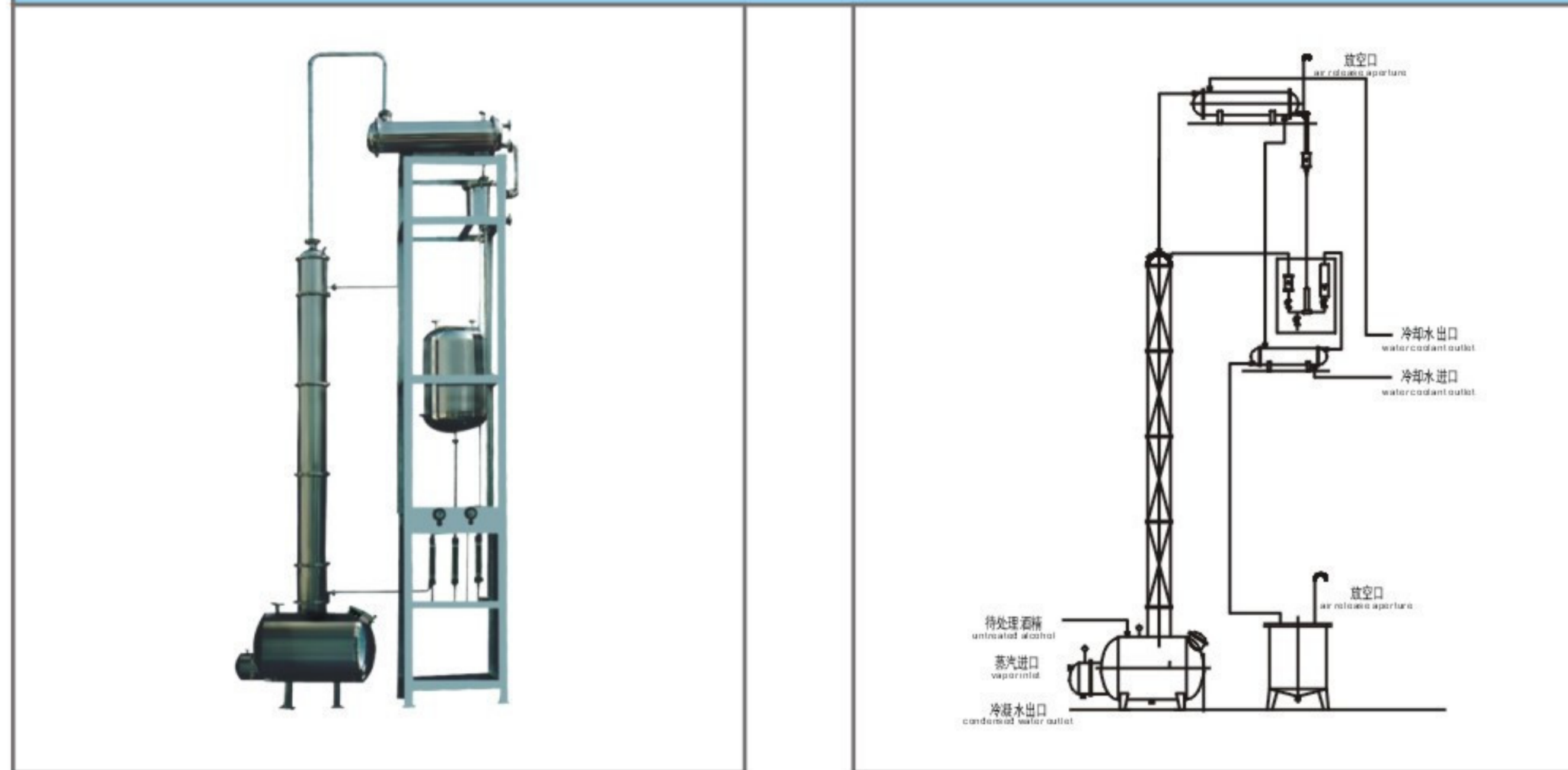
JH-series Alcohol Recovery Tower is made up of tower pot, tower body, condenser, cooler, and high-level storage tank, which is applicable to pharmacy, food, light industry and chemical engineering works for thin alcohol recovery. The parts of this equipment that contact materials are all made of stainless steel SUS304 or SUS316L, so the equipment has good corrosion-resistant capability and can save energy, protect environment, reduce production cost and increase efficiency. This device can distill thin alcohol of 30~50 degree to 93~95 degree. The residual liquid has low discharge alcohol content, which is in accordance with environment protection requirement. The device adopts highly-effective stainless steel corrugated loading material. The parts contacting material are all made of stainless steel. The equipment surface has got surface polishing treatment. GMP medical standards.

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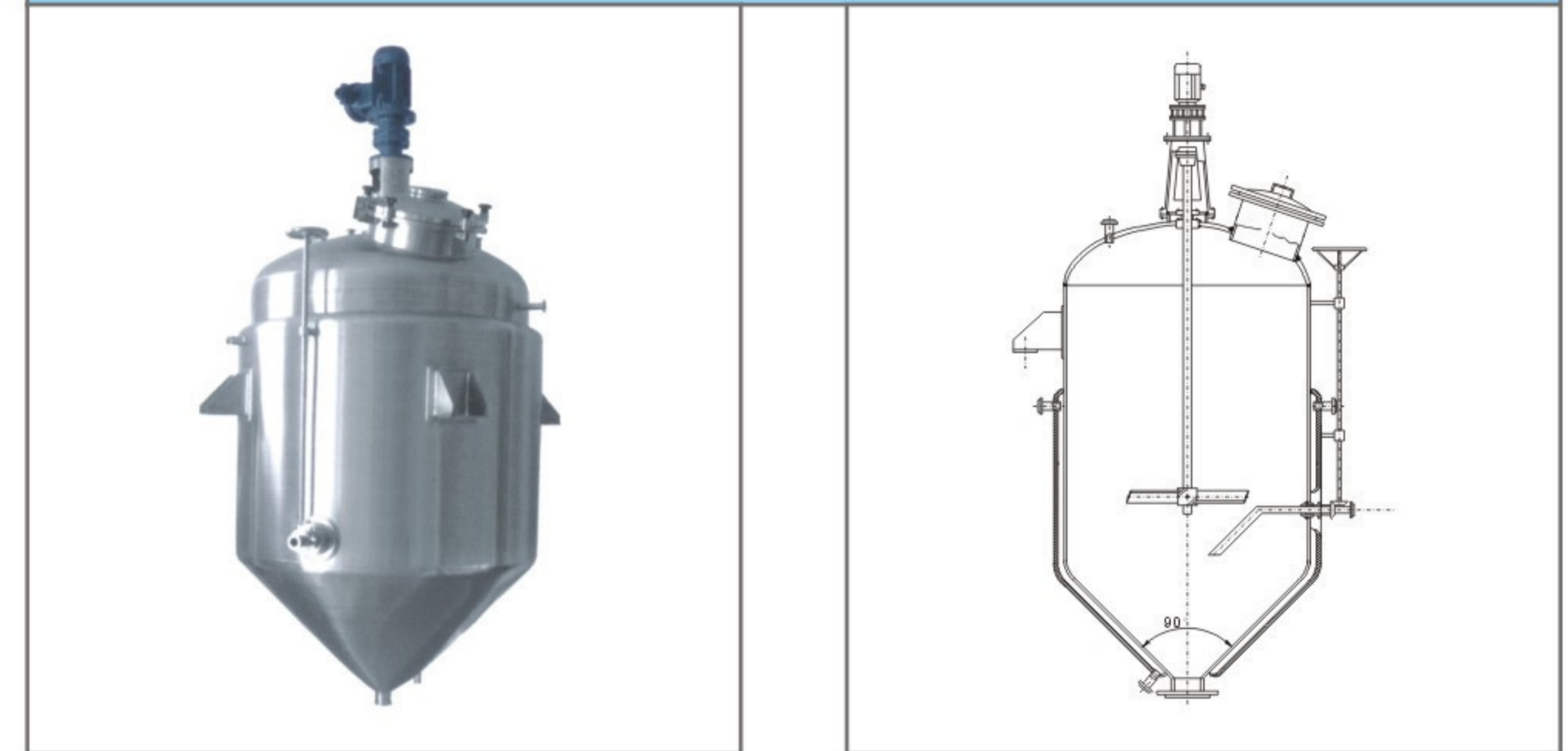The device is suitable for alcohol deposition in such industries as China traditional drug, oral liquid, foodstuff, health product, chemistry. The alcohol deposition tank is a special device for alcohol deposition, also titled solvent change or alcohol bar. It is mainly used to freeze the solution from concentration of water-decocted China traditional drug or make alcohol deposition under normal temperature, also to execute the water deposition of the concentrated liquid after the alcohol deposition of China traditional drug. The device includes the ellipse seal head with clip sleeve, tapered-shaped base of cylinder, ommer-installed three-paddle miller and special finely-regulated revolve-out liquid pipe, etc. The mode of water-decocting and concentration of China traditional drug as follow: First fill the alcohol to form mixture, secondly under low temperature freeze to make deposition to separate the solid from the liquid to improve the purity and cleanness of the extraction liquid of China traditional drug. The device is made with stainless steel of SUS304. The high pressure automatic spray & cleaning system is available. It is available to fill the low temperature of cooling water through the clip sleeve. The mechanical seal in the transmission part and the explosion-proof motor can ensure the security of the production process, meeting with the GMP standard. The surface is treated with polish or surface polishing.

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Usage

The equipment is applicable for the filtration operation for such industries as traditional Chinese medicine, food and chemical industry etc.

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Working principle

The filtration method is to add continuously leaching solvent into the rough powder of medicine material to make it filtrate the medicine powder, and then the leaching liquid will flow from the bottom outlet. During the filtration, the solvent will filter into the cell of medicine material to dissolve large amount of soluble matter. The concentration increases, and it will move downward after the density increases. The leaching solvent at the upper layer or the thin leaching liquid will exchange the position and produce good concentration difference to make the scattering be carried out smoothly. Thus, the leaching effect is better than the dipping method, and the extraction will be comparatively complete.

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The equipment is composed of tank body, taper body, oval seal head (or flat cover), pneumatic slagging door and pneumatic operation bench etc. It has such characteristics as equal solvent distribution and convenient slagging etc.

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PYG series processing Tank (Reactor)

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The liquid arrangement tank strictly following with the 31st and 32nd clauses of 《Medicine Production Quality Management Norm》-GMP General Principles (1998 Revised Edition), and is designed, manufactured and accepted according to the requirements of GB150-98 《Steel Pressure Vessel》 and JB/T4735-97 《Steel Welding Normal Vessel》.

- ◆ Isolation material of the vessel adopts high-temperature resistant rock wool and anti-aging PU foam.
- ◆ ISO clamp, inner container adopts imported stainless steel SUS316L or SUS304, internal surface is polished to $Ra \leq 0.28 \mu m$, and external surface polishing mirror polish, sand blasting or cold-rolled nature-color surface polishing. All is in accordance with GMP medical standards.
- ◆ Sealing of mixer adopts imported sanitation-grade mechanical stirring enclosing or magnetic totally-enclosed stirring device. The mixing rotary speed is optional.
- ◆ Orifice is set with level gauge (static-pressure type, capacitance type, non-contact type, ultrasonic type and glass tube type), air respirator, thermometer (digital display type or dial plate type), CIP cleaning mouth, viewing mirror, explosion-proof viewing lamp (viewing mirror and viewing lamp integration), SIP sterilizing mouth, liquid entrance-exit gate and sanitary inlet.
- ◆ Volume Specification: 100L-10000L or to be designed and manufactured at customers' actual request.
- ◆ Relevant GMP testing data are attached (including material report, external components certificate of quality and verification form).

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
FJ insulation and Cooling Fermentation Tank (Seed tank, crystallizing tank)

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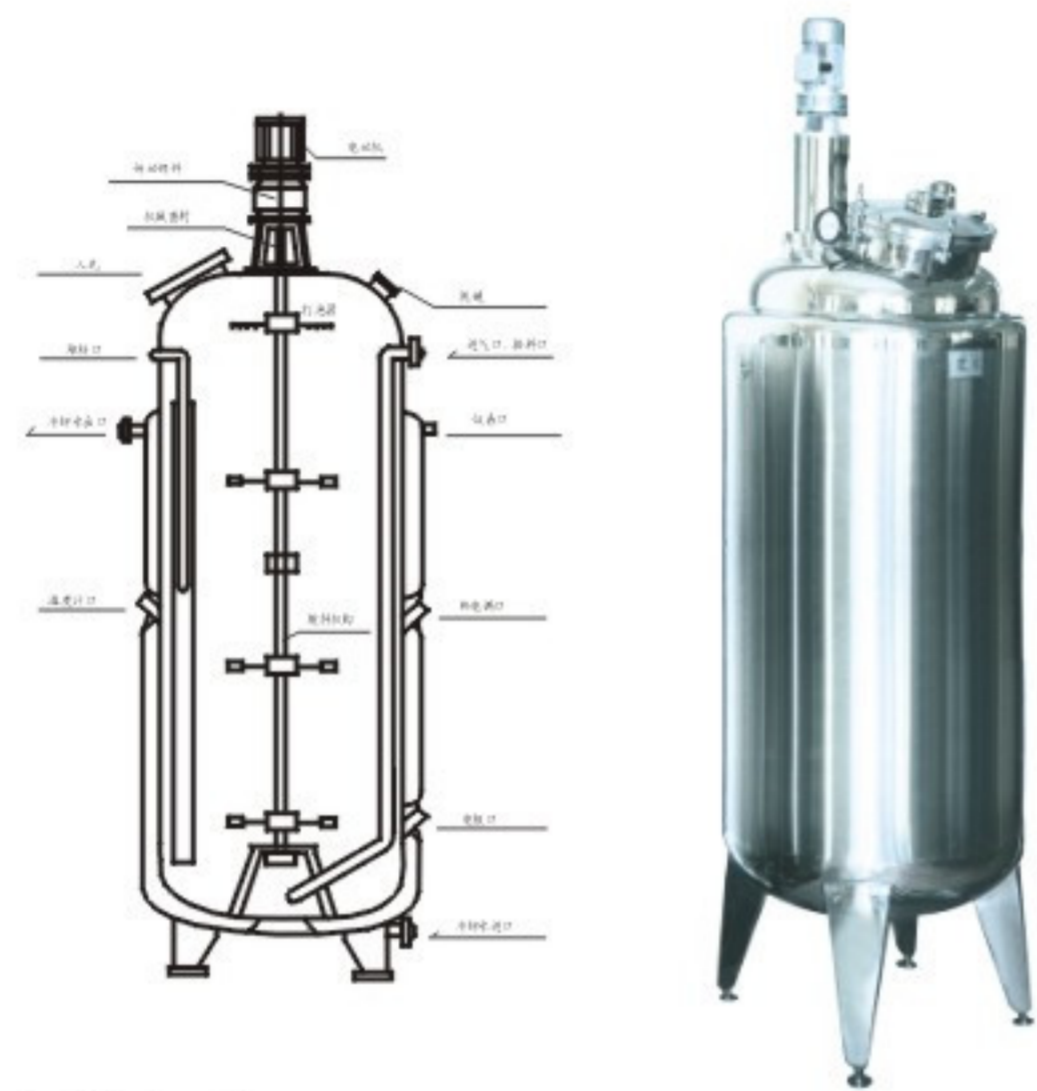Fermentation Tank(seed tank,crystallizing tank) is widely applied in pharmacy, fine chemical engineering and biological engineering in dustries. The tank body has interlayer and heat-insulating layer (for heating, cooling or heat insulating) The tank body , upper and lower closure heads(or taper) all adopt spinning R-angle processing. The internal wall of the tank has got mirror polishing with finish Ra≤0.4 μm. It has no sanitary dead corner and its wholly-enclosed design ensures that materials are always. Mixed and ferment at non-pollution state. It is equipped with devices including air respirator, cleaning ball, sanitation inlet , viewing mirror, and viewing lamp. The fermentation process of Pharmacy, fine chemical engineering and biological products is aseptic and untainted. Fermentation tank adopts aseptic system which helps avoid microorganism pollution in the air and greatly extends quality warranty period and purity of the products. We specially designed and installed aseptic respiratory air hole or aseptic normal-pressure fermentation system, which can finish circulating heat-up and cooling through medium heat-up and cooling. The parts that contact materials are all made of stainless steel SUS304 or SUS316L. Heat preservation adopts polyurethane blister or pearl cotton. External surface adopts sand grinding surface polishing treatment. All is in accordance with GMP standard Volume : 100-20000L.

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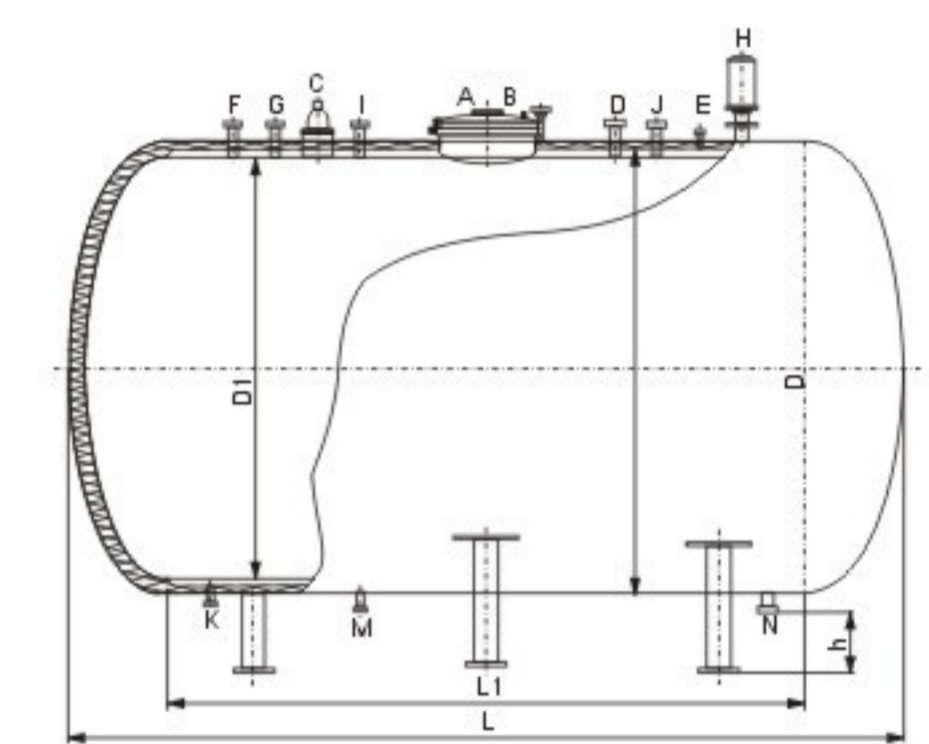
- ◆Type: horizontal type or vertical type
The horizontal distilled water tank series strictly follow with the 31st and 32nd clauses of 'Medicine Production Quality Management Norm--GMP General Principles' (1998 Revised Edition), and is designed, manufactured and accepted according to the requirements of Gb150-98 'Steel Pressure Vessel' and JB/T4735-97 'Steel Welding Normal Vessel'.
- ◆Isolation material adopts high-temperature resistant rock wool or anti-aging PU foam.
- ◆ISO clamp, inner container adopts SUS316L or SUS304, internal surface polished to Ra≤0.28 μm, and external surface polishing, mirror polish, sand blasting or cold-rolled nature-color surface polishing. All is in accordance with GMP.
- ◆level gauge (static-pressure type, capacitance type, non-contact type, ultrasonic type and glass tube type), air respirator, thermometer (digital display type or dial plate type), CIP cleaning mouth, SIP sterilizing mouth, liquid entrance-exit gate and sanitary inlet.
- ◆Volume Specification: 100L-30000L or to be designed and manufactured at customers' actual request.
- ◆Relevant GMP testing data are attached (including material report, external components certificate of quality and verification form).

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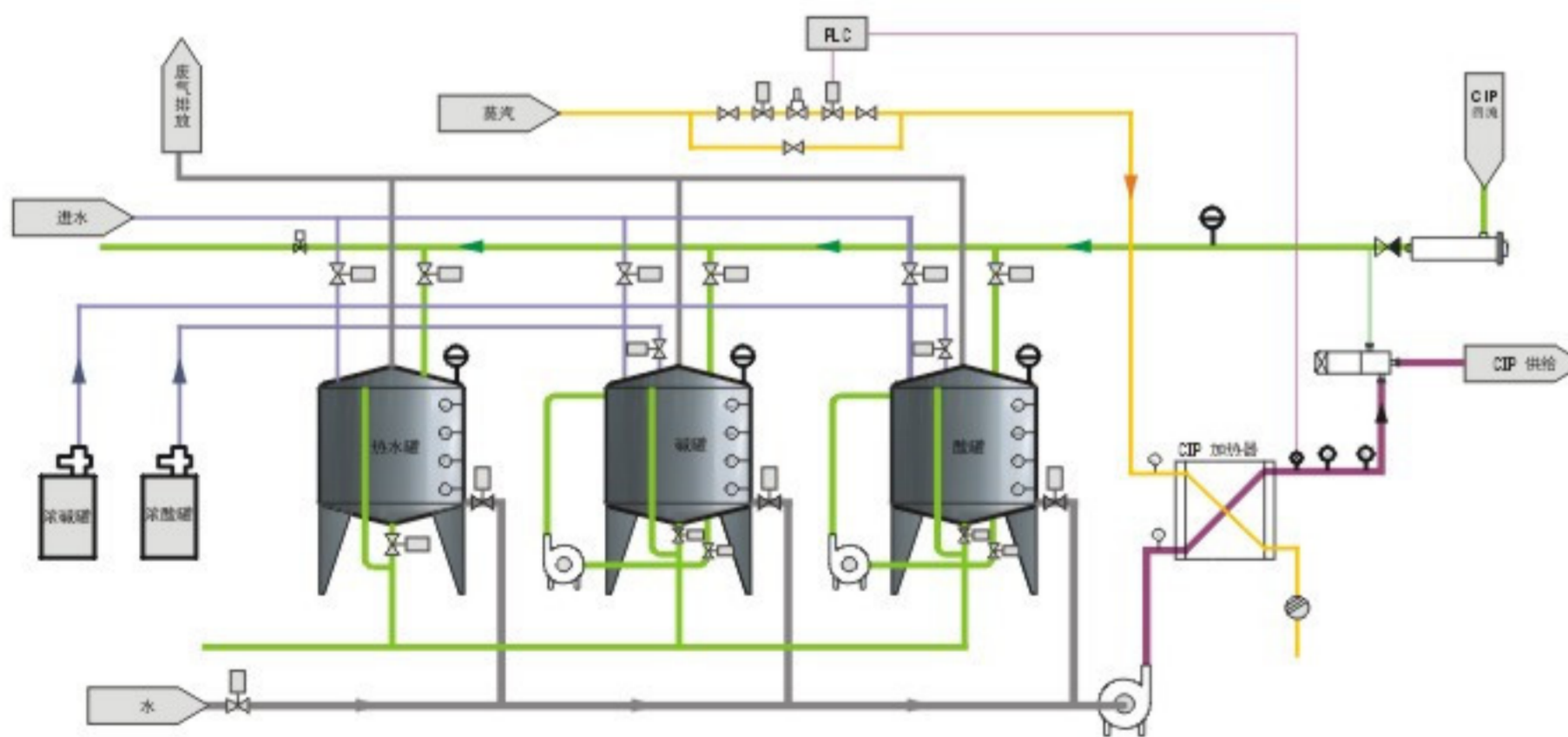
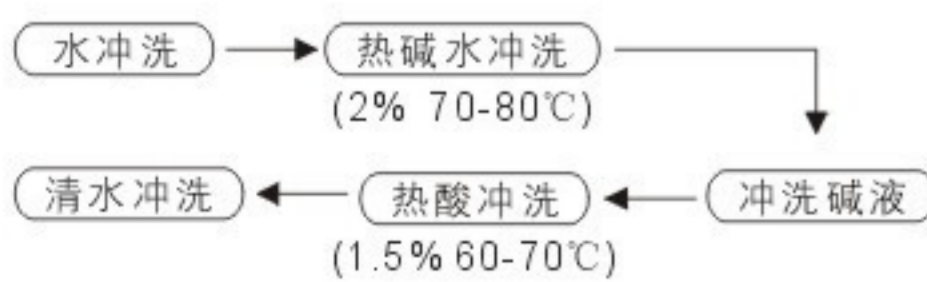
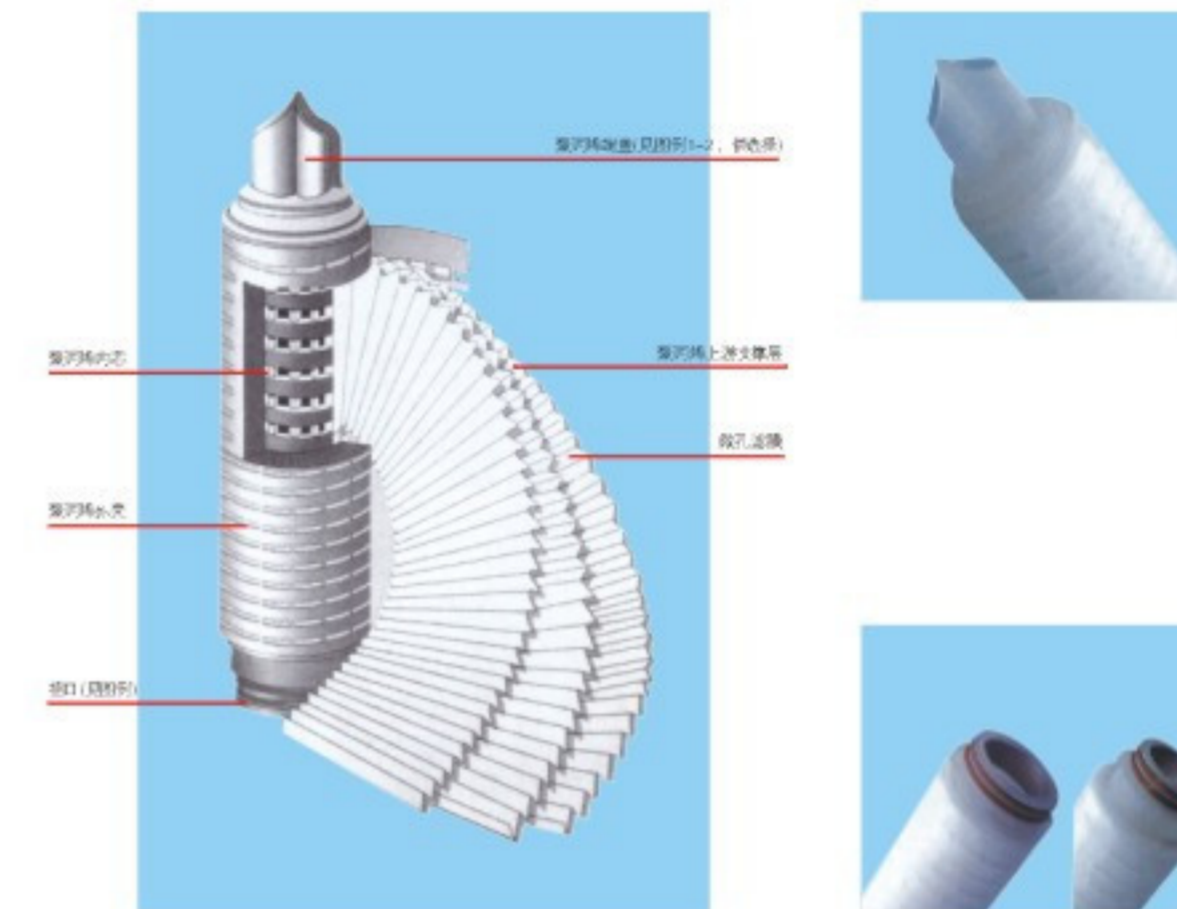
Microporous membrane barrel filter

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- ◆Material: SUS304 or SUS316L
- ◆Application: Winery and brewery, beverage, juice, vinegar, pharmaceutical, chemistry, etc
- ◆filter cartridge Qty: 1pcs, 3pcs, 5pcs, 7pcs, 9pcs, 11pcs, 13pcs, 15pcs
- ◆size: 10", 20", 30", 40"
- ◆Cartridge material: EPDM, PP, nylon, etc
- ◆Accuracy: 0.1 μm, 0.22 μm, 0.45 μm, 1 μm, 3 μm, 5 μm, 10 μm,

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- ◆Pump head material: SUS304 or SUS316L
- ◆Application: For dairy, food, Juice, pharmaceutical and biochemistry industry delivery.
- ◆type : Portable or Fixed
- ◆impeller: Close type of open type
- ◆Motor: Option (Chinese type or European type)
- ◆Mechanical seal are not connect with material
- ◆pump heat drainage hygiene design ,no dead angle
- ◆Max lift: Option in the reasonable range

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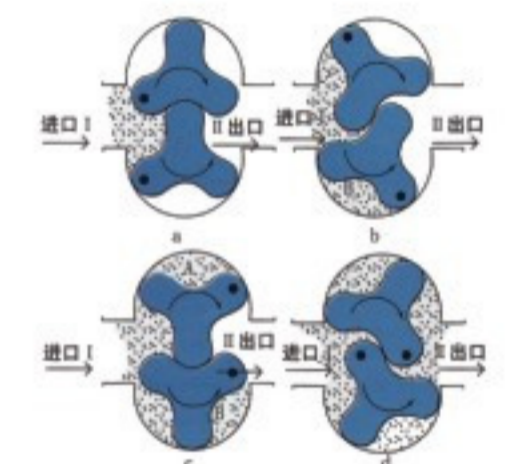
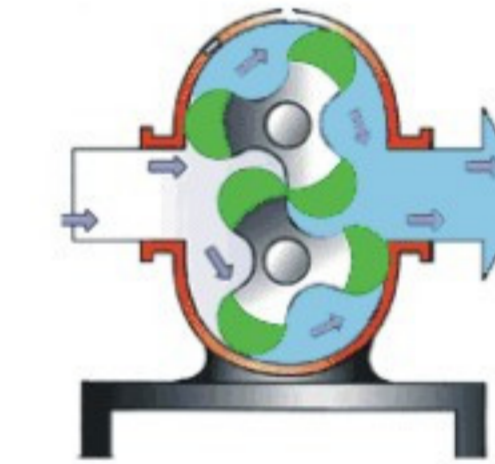
- ◆Capacity: 500L/h ~14,000 L/h
- ◆Pump head material: SUS304 or SUS316L
- ◆Application: For dairy, food, Juice, pharmaceutical, cosmetics, detergents, hair gel, aromatic oils and biochemistry industry delivery.
- ◆type : (Portable or Fixed)
- ◆pump head: Horizontal or vertical delivery
- ◆Motor: Option (Chinese or European type) Variable speed, variable frequency, fixed speed gear
- ◆configure: Mechanical seal, O ring
- ◆rotor: unifoliate, bifoliate, trifoliate
- ◆material temp range: -30°C~ 150°C
- ◆other element: Processing liquid viscosity

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双桶过滤器
Double-barrel Filter

列管式浓缩器
Pipe Concentrator

盘管式真空浓缩锅
Tubular vacuum concentration pot

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Sanitary stainless tank system configuration

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The pipe and valve are highly effect the hygeine demand of the equipments, which are one of the most important demand of the equipments.so we make each detail of the fittings, material are strictly following SUS304 SUS316L , polished $Ra \leq 0.28 \mu m$, outside finished as matt polish, we have fitting in different standard: ISO, DIN, IDF, SMS, RJT etc, GMP standard.